



BIO-HAZARD SPILL KIT 5 Ltr

SPLSK05BHZ ABSORBENCY : 05 Ltr

CONTAINS :

- › One Set of Personal Protective Equipments
 - › Safety Goggles : 1 Nos
 - › 3Ply MB Mask : 1 Nos
 - › Safety Gloves : 1 pair
- › Clinical Wastage Bag with Tie : 3 Nos
- › Absorbent Granules : 500 Gms
- › Caution Mat Yellow : 2 nos
- › Chemical Absorbent Pads (Green & Blue) : 6Nos
- › Chemical Absorbent Socks (8cm x 1.2m) : 1 Nos
- › MAT Bag 15 Ltr Capacity : 1 Nos

SPILFIX SPLSK05BHZ Biohazard Spill Kits Specially Customized design and packed for quick & Efficient Response against Small Spillage.

- › Put on Protective Gears Supplied with Kit.
- › Sprinkle Granular over the Spill and leave for 2 min.
- › Once Granules have absorbed the spill, collect same using x-ray sheet & Scoop and brush.
- › Sanitise the Spill area with disinfectant liquids.
- › Clean the Spill area using Mats & Pads.
- › Bag and contents must be disposed of in accordance with local waste disposal policies for biohazard waste.
- › As per NABH standard and satisfied ISO guidelines