



BIO-HAZARD SPILL KIT 25 Ltr

SPLSK25BHZ ABSORBENCY : 25 Ltr

CONTAINS :

- › One Set of Personal Protective Equipments
 - › Safety Goggles : 1 Nos
 - › 3Ply MB Mask : 1 Nos
 - › Chemical Coverall : 1 Nos
 - › Safety Gloves : 1 pair
 - › Baricadding Tape : 1 Nos
 - › Shoe Cover : 1 Pair
- › Clinical Wastage Bag with Tie : 5 Nos
- › Absorbent Granules : 2500 Gms
- › Caution Mat Yellow : 3 nos
- › Chemical Absorbent Pads (Green & Blue) : 20Nos
- › Chemical Absorbent Socks (8cm x 1.2m) : 2 Nos
- › Pan & Brush : 1 Nos
- › SS 304 BOX WITH LUCK : 1 Nos
- › Rejected X-Ray Sheet : 2 Nos
- › Disinfectant Liquids : 250 ML
- › Blotting Papers : 5 nos

SPILFIX SPLSK100BHZ Biohazard Spill Kits Specially Customized design and packed for quick & Efficient Response against Large Spillage.

- › Put on Protective Gears Supplied with Kit.
- › Sprinkle Granular over the Spill and leave for 2 min.
- › Once Granules have absorbed the spill, collect same using x-ray sheet & Scoop and brush.
- › Sanitise the Spill area with disinfectant liquids.
- › Clean the Spill area using Mats & Pads.
- › Bag and contents must be disposed of in accordance with local waste disposal policies for biohazard waste.
- › As per NABH standard and satisfied ISO guidelines